

KARL FISCHER TITRATOR WITH 21 CFR PART 11 COMPLIANCE

KAFI SMART



LABINDIA
ANALYTICAL
WHERE COMMITMENT IS CULTURE

Automatic Karl Fischer Titrator & KAFI⁺

Karl Fischer titration is an accurate, rapid and efficient method for determining the water content in samples. When the sample is titrated in the presence of SO_2 , I_2 , and an organic base, moisture from the sample extracted in the solvent, can be quantitatively estimated. Dry methanol is used as a solvent in Karl Fischer titration.

The indicator-electrode is a two-pin platinum electrode across which a constant current is generated by a polarizing current source. For starting, very low volume of methanol as a solvent is needed in the specially designed vessel.

Background signal correction is done through an offset mode.

Calculated doses of Karl Fischer reagent are added to the solution in a hermetically sealed vessel as per stop criteria selected and the end-point volume is evaluated automatically. The instrument continuously monitors moisture leak in the vessel and neutralizes it automatically.

The built-in reagent standardization procedure by water (H_2O) or Sodium Tartrate ($\text{C}_4\text{H}_4\text{Na}_2\text{O}_6 \cdot 2\text{H}_2\text{O}$) for titre factor/concentration (F) is incorporated.

The liquid handling path comprises teflon tubing, teflon-lined valve, and a gastight syringe with teflon plunger. It creates a chemically inert system for this highly sensitive analysis.

The magnetic, adjustable-speed stirrer is a part of the system in which the sample is stirred vigorously to extract the moisture efficiently.

The result, including the leak rate, is printed in a tabulated form and is also displayed on back lighted liquid crystal display (LCD) screen.

MODES OF OPERATION

Karl Fischer titration is a dead stop titration process which monitors the mV drift to be lower than the stop band, and the end of the titration is detected when the leak rate is close to or less than stop band. The leak rate is displayed and also printed. There are two main modes of operation:

- (a) In minimum dose mode where the end dose is same as the minimum dose, the titration will be stopped when the mV does not rise above the stop band in specified delay time.
- (b) In drift mode, the titration will be stopped if the calculated drift value is less than or equal to entered drift value in $\mu\text{l}/\text{min}$.

APPLICATIONS

- Pharmaceutical
- Food and Beverages
- Petrochemicals
- Cosmetics
- Organic / Inorganic Chemicals etc.

FEATURES


- Advanced Microcontroller based user-friendly, state-of-the-art product design with User interactive software for ease of operation with protection against invalid entries.
- Quick interchangeable imported burette assemblies with intelligent recognition for its volume size. Burette validation factor for dispensing correction is available for true end point volume.
- System recognizes proper connectivity of other peripherals like Burette, Stirrer, Electrode, Pen Drive etc. Gives indication incase of improper connectivity.
- Large memory capacity for method storage with suitable scientific parameters having GLP compliance.
- Sample Name & Identification Number with Date and Time for authentication. Daily Auto Incremented Run number and Factory entered CUSTOMER NAME & Instrument Sr. No. on report printouts make the system foolproof and GLP compliant.
- Quick monitoring, and automatic neutralization of moisture leak into vessel to keep it ready for next titration.
- End point delay up to 100 sec for slow moisture releasing samples.
- On line leak rate correction available.
- Microcontroller based variable speed, magnetic stirrer with digital indication.
- Calculation modes :
Weight /Weight
Volume/Volume
Volume/Density
- Selectable report format, complying with GLP requirements :
A) Report giving titration parameter and result.
B) Report of method parameters.
C) Condensed report of titration parameter and result.
D) Graph printout of reaction trend

Reports can be obtained even after resetting / power off / power failure conditions.

- Statistic function with run selectivity for finding Mean, S.D., R.S.D. and C.V. of last 10 repeat run results could be viewed or printed.
- ASTM standard compliant for analysis of oil samples.
- Date & Time display and report printout with run time indication.
- Balance interface to directly transfer the sample weight.

21 CFR PART 11 COMPLIANCE

- Audit Trail for all activities with search facility, report generation and printing
- 200 User ID's with alphanumeric entries of user name, password and role based privileges selection
- Multi-level roles with password protection
- User authentication is performed for each and every operation done by user
- Customizable PDF report file can be created through print
- USB Printing eliminates the need of serial port to connect with instrument
The user can take printout on printer via WiFi connectivity
- Electronic signature functionality
- Manual Archive and Data Backup facility available

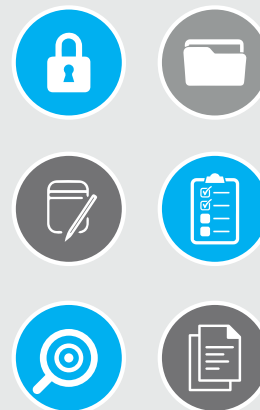

Optional

- IQ, OQ, PQ, documents available
- Auto Fill & Drain Pump (i.e. for draining out waste from reaction vessel & to fill with fresh solvent)

SPECIFICATIONS

- **Principle** : Karl Fischer method of volumetric water determination.
- **mV range** : ± 3200 mV.
- **Accuracy** : ± 1 mV.
- **Polarized Measurement range** : 0 to ± 3200 mV
- **Polarized sensor Resolution** : 0.1 mV
- **End point detection** : Voltametric.
- **Polarising Current Range** : 1 μ A to 80 μ A in 8 steps.
- **Polarization Current Resolution** : 0.1 μ A
- **Polarization Current Accuracy**: 1.0 μ A
- **Cut-off criteria** : Delay or drift - user selectable.
- **Measuring Range** : 10 μ g to 500 mg (10 ppm - 100%)
- **Measuring Accuracy**: Better than 0.3% of respective Volume
- **Repeatability** : Better than 0.3% at > 10 mg H₂O
(at ideal analytical conditions)
- **Sensor** : Dual pin platinum electrode.
- **Sensor input** : Through special TNC connector.
- **Burette** : 5ml or 10ml capacity, interchangeable, with auto recognition.
- **Burette Resolution** : 1/5000 for 5 ml (i.e. 1 microlitre),
1/10000 for 10 ml (i.e. 1 microlitre)
- **Filling time** : < 20 sec.
- **Stirrer** : Magnetic capsule type, microcontroller based speed control with digital indication.
- **Display** : 7" TOUCH SCREEN
- **Method Storage** : 16 GB data storage capacity for 200 methods & results.
- **Results** : mg/ml, % , ppm, mg/g and on-line leak rate with cumulative titrant consumption.
- **Drift Measurement**: Online, $< 2\mu$ g/min

REPORT FORMAT :

- a) **Method parameter.**
 - b) **Titration analysis report.**
 - c) **Titration analysis condensed report.**
 - d) **Statistics and on-line leak rate report.**
 - e) **Graph for reaction trend**
-
- **Input/Output peripheral interface** :
 - a) **USB Port** : 1 No. for printer / pen drive
 - b) **Serial Port** : 1 No. for Balance.
 - **Power requirement** : 230 V AC ± 10 % , 50 Hz.
 - **Environmental Operating Conditions** :
 - a) **Operation** : Indoor
 - b) **Temperature** : Ambient to 45 °C
 - c) **Humidity** : 5 to 90% non-condensing.

LABINDIA

A N A L Y T I C A L

WHERE COMMITMENT IS CULTURE

www.labindia-analytical.com / For Enquiries: sales.mfd@labindia.com

Labindia Analytical Instruments Pvt. Ltd.

HEAD OFFICE & REGIONAL OFFICES:

Thane: 201, Nand Chambers, LBS Marg, Near Vandana Cinema, Thane (W) - 400 602.
Tel.: +91-22-2598 6000 / 2598 6110 Fax: +91-22-2541 0420 / 2533 5940
E-mail: sales.mfd@labindia.com

Delhi: G-4, Pal Mohan Sadan, 26/32, East Patel Nagar, New Delhi - 110 008
Tel.: +91-11-4330 6001 / 10 Fax: +91-11-2585 1066

Chennai: B-1, Alsa Regency, 165, Eldams Road, Chennai - 600 018.
Tel.: +91-44-2434 7008 / 2432 0352 Fax: +91-44-2434 6328

Kolkatta: 165-A, S. P. Mukherjee Road, Kolkatta - 700 026.
Tel.: +91-33-2466 3362 Fax: +91-11-2466 1352

BRANCH OFFICES:

Bangalore: #105, Bhimajyothi Colony, West of Chord Road, Basaveshware Nagar,
Above Coffee Day, Bangalore - 560 079 Tel.: +91-80-2323 0919 / 20 / 21
Fax: +91-80-2323 0924

Lucknow: 403, 4th Floor, Sahara Shopping Complex, Faizabad Road, Indira Nagar,
Lucknow - 226 016. Tel.: +91-522-2346 535 / 2346 496 Fax: +91-522-23488 847

Hyderabad: 6-3-1090/1/, Uma Hyderabad House, 2nd Floor, Somajiguda, Raj Bhavan
Road, Hyderabad - 500 082. Tel.: +91-40-4444 3456
Fax: +91-40-2331 3312

Tiruvanthapuram: Megha Complex, Ground Floor, TC No. 26/857, Women College Road,
Thycaud, Thruvanthapuram - 695 014 Tel.: +91-471-2324 064 / 2320 082
Fax: +91-471-2320 082

Vadodara: 605, Imperial Heights, Akshar Chowk, Old Padra Road, Vadodara - 390020.
Tel.: +91-265-2986 005 / 2986 006

Pune: Office No. 25, 3rd Floor, Lokmanya House, near CNG Petrol Pump, Paud Road,
Kothrud, Pune - 411 038 Tel.: +91-20-2538 0090

Chandigarh: S.C.O. 208-209, 4th Floor, Basera Building, Sector-34 A,
Chandigarh - 160 022, U.T. Tel.: +91-172-4090 001-07 / 4090 009
Fax: +91-172-4090 008

FACTORY:

Plot No. EL-72, Electronic Zone, TTC Industrial Area, Thane Belapur Road,
Navi Mumbai - 400 705.
Tel.: +91-22-2762 6660 / 61 / 62 / 63 / 6516 3480 / 81 / 6107 6666
Fax: +91-22-27626664 E-mail: lipifactory@vsnl.net

APPLICATION & TRAINING CENTRE:

R-909, TTC Industrial Area, Thane Belapur Road, Rabale, Navi Mumbai - 400 701.
Tel.: +91-22-2760 6955 / 6107 6555 Fax: +91-22-2760 6706



Striving to become the best individuals, we endeavour to foster the best team. Performing sensibly, we try to achieve the best efficiency. Working innovatively, we seek to make the best products. Listening patiently, we excel to offer the best service. So, no matter what you needs are, **come to us, GET THE BEST**

LABINDIA reserve the right to change specification without notice as part of its continuous programme of product development.